





Scherzo SW-C18

100 x 3 mm

Application

LC-MS: Antipsychotics

抗精神薬のLC-MS分析

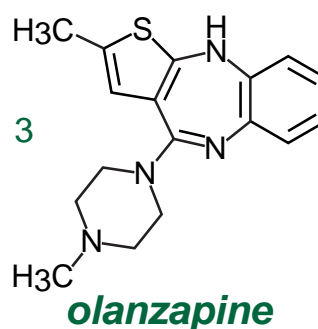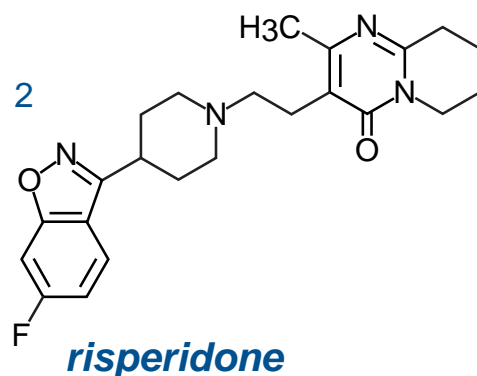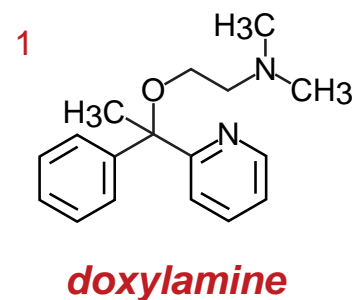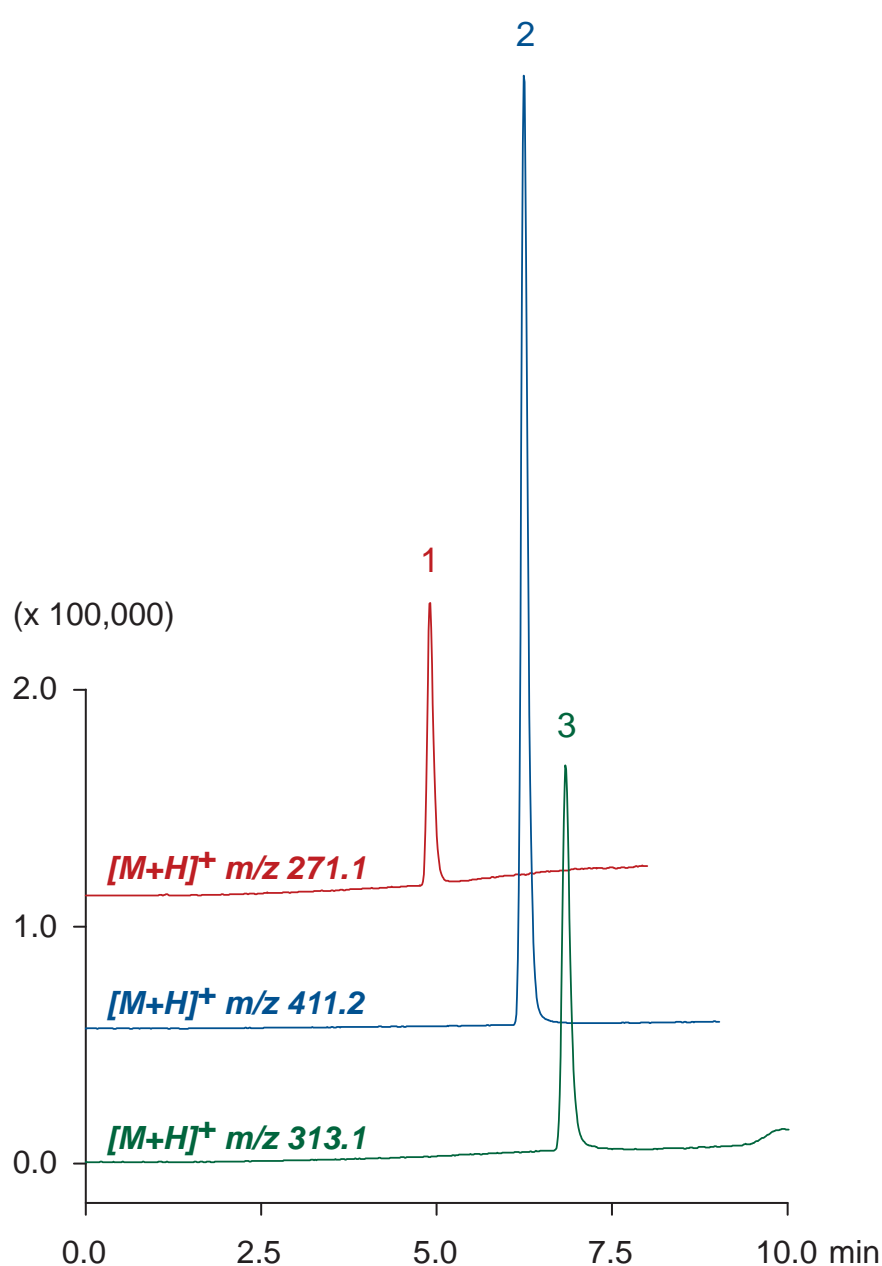

Scherzo SW-C18, 100 x 3 mm
 A: 10mM ammonium formate
 B: acetonitrile / 50mM ammonium formate = 50 / 50
 50-100%B (0-10 min), 100%B (10-15 min)
 0.4 mL/min (6 MPa), 37°C,
 1 µL (1µg/mL, 10mM ammonium formate)
 Single Quad. MS (ESI)

Distributor

LCC Engineering & Trading GmbH, Steinbruchstr. 4, CH-4622 Egerkingen, Switzerland.
 www.chromatographyshop.com , info@chromatographyshop.com,
 Tel: +41 62 398 52 71 , Fax: +41 62 398 52 74