




Scherzo SW-C18
Scherzo SS-C18

50 x 3 mm

Application

LC-MS: Glutamyl-Valyl-Glycine

グルタミンバリルグリシン(コク味ペプチド)のLC-MS分析

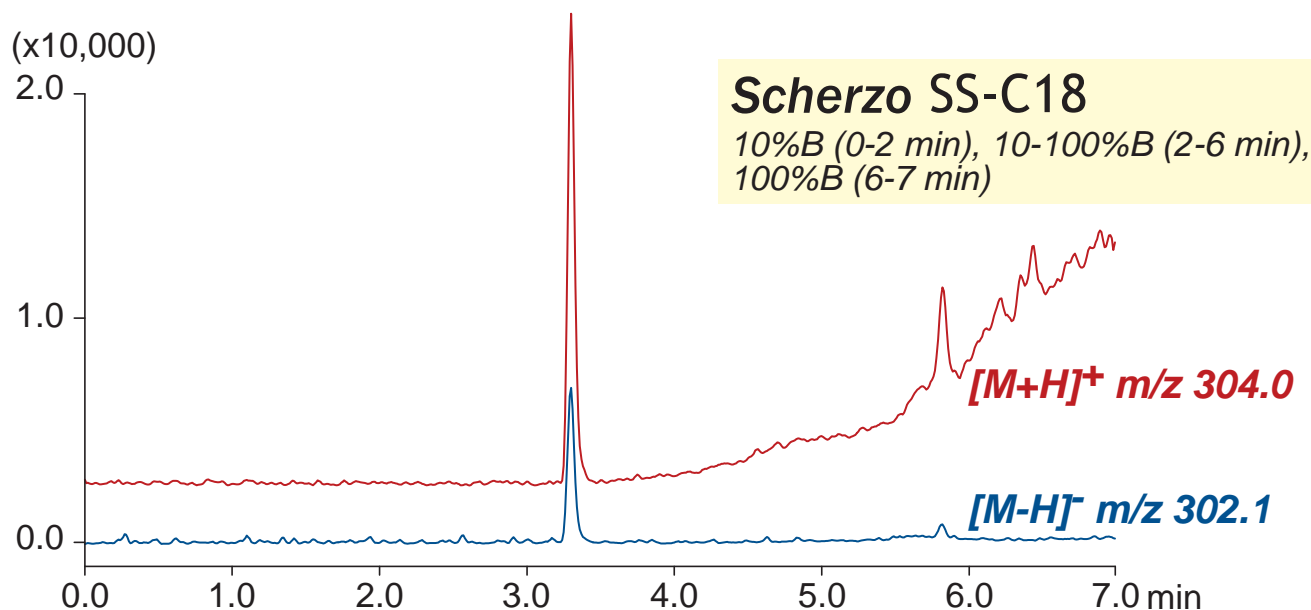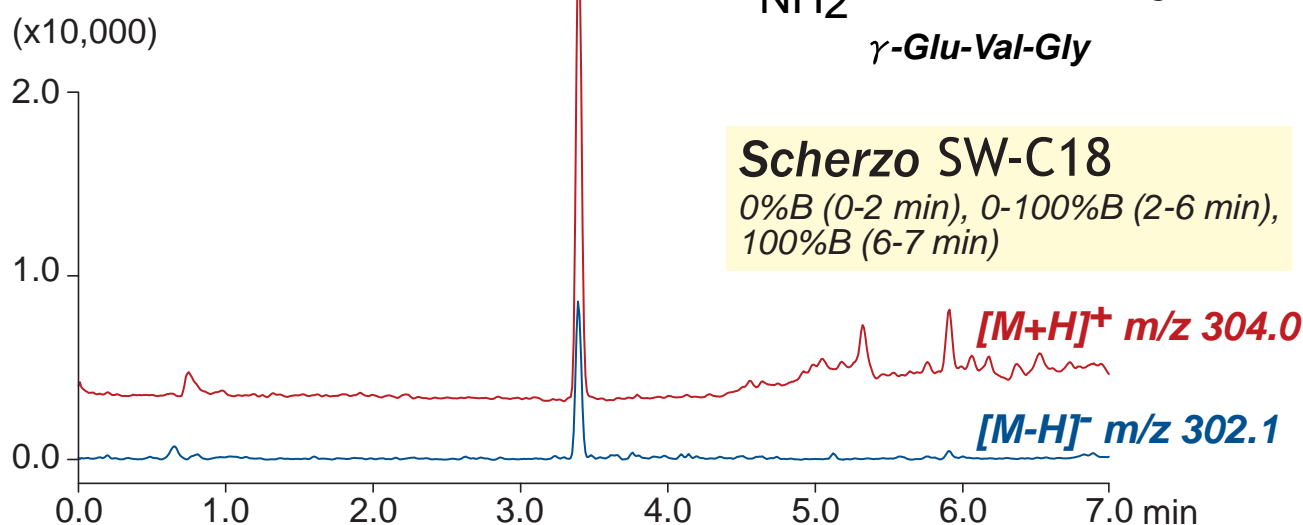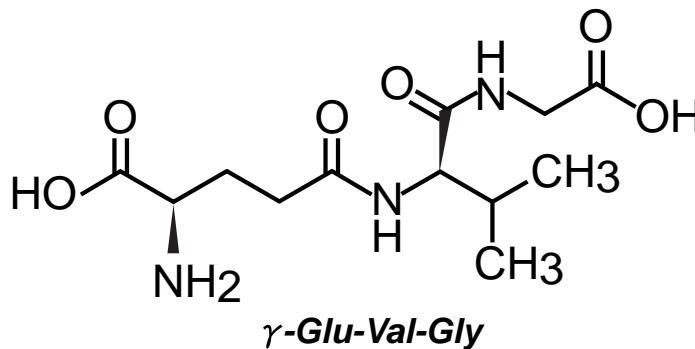

50 x 3 mm

A: water / formic acid = 100 / 0.1

B: acetonitrile / 100mM ammonium formate = 40 / 60

0.4 mL/min (4 MPa), 37°C, 20 μ L (500ng/mL, water / HCOOH = 100 / 1)

Single Quad. MS (ESI)

Distributor

LCC Engineering & Trading GmbH, Steinbruchstr. 4, CH-4622 Egerkingen, Switzerland.

www.chromatographyshop.com , info@chromatographyshop.com,

Tel: +41 62 398 52 71 , Fax. +41 62 398 52 74