


Conditions

Column: ACE 3 AQ
Dimensions: 150 x 3.0 mm
Part Number: ACE-116-1503
Mobile Phase: A: 0.1% formic acid in H₂O
 B: 0.1% formic acid in MeCN
Gradient:

Time (mins)	%B
0	0
10	100

Flow Rate: 0.4 mL/min
Injection: 5 µL
Temperature: 30 °C
Detection: Exacte Orbitrap high resolution MS
 ESI positive ion mode
 Capillary temperature: 350 °C

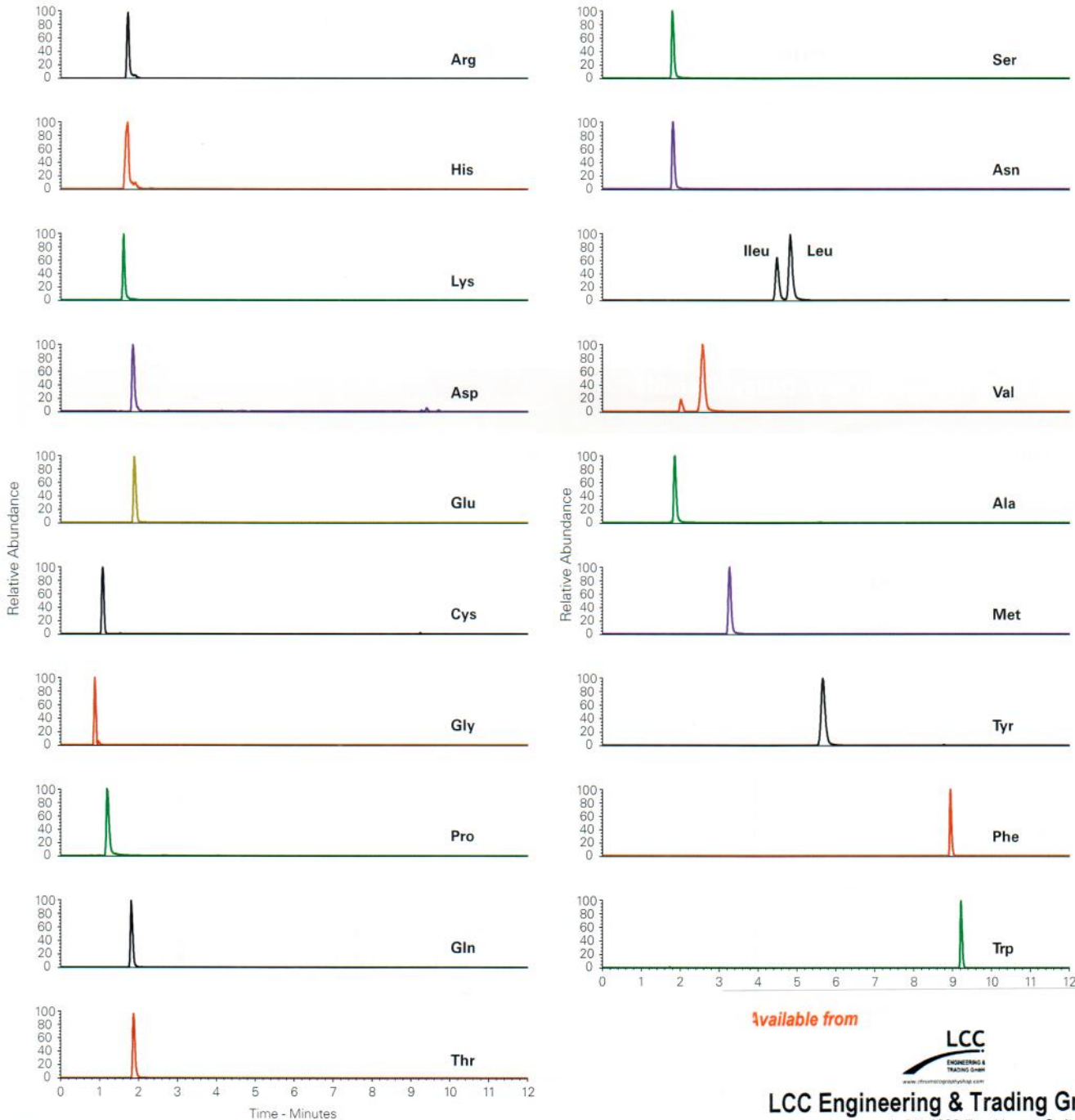

Available from



LCC Engineering & Trading GmbH
 Steinbruchstrasse 4, CH-4622 Egerkingen, Switzerland
 Tel.: +41 62 398 5271, Fax.: +41 62 398 5274
info@chemsupply.ch
www.chromatographyshop.com

Reproduced with permission of Fera Science Ltd, York, UK