


Scherzo SM-C18

50 x 3 mm

Application

LC-MS: biogenic polyamines

生体ポリアミンのLC-MS分析

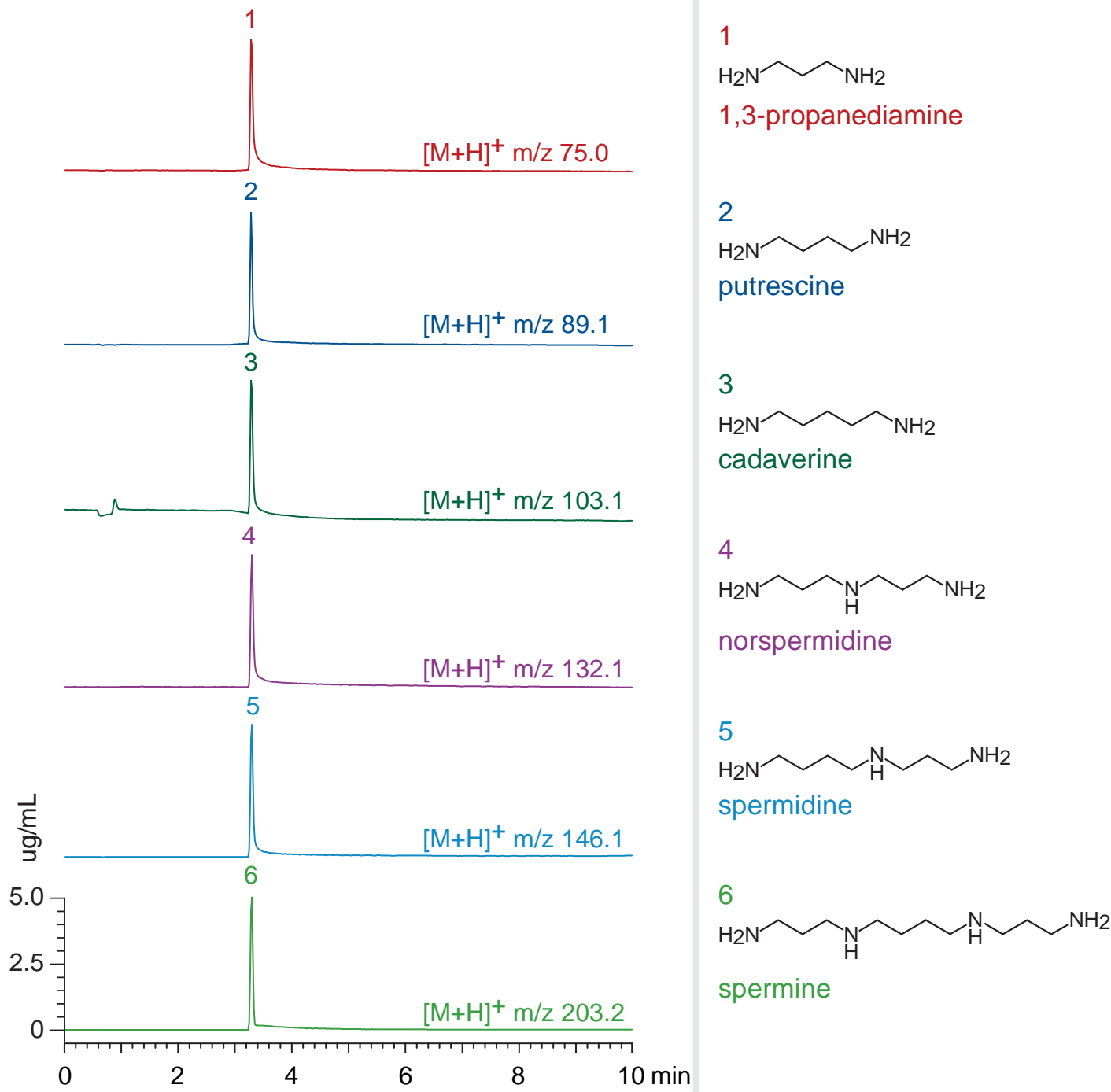

Scherzo SM-C18, 50 x 3 mm

A: 5 mM ammonium formate

B: acetonitrile / water / formic acid = 20 / 80 / 0.8

0 %B (0-1min), 0-70 %B (1-8min), 100 %B (8-10min)

0.4 mL/min (4 MPa), 37°C, 10uL (5ug/mL, 50mM AcONH<sub>4</sub>)

Single Quad. MS (ESI, positive)

**Distributor**

LCC Engineering & Trading GmbH, Steinbruchstr. 4, CH-4622 Egerkingen, Switzerland.

www.chromatographyshop.com, info@chromatographyshop.com,

Tel: +41 62 398 52 71, Fax: +41 62 398 52 74