



Scherzo SM-C18

100 x 3 mm

Application

LC-MS: biogenic amines and precursor amino acids

腐敗アミンと前駆アミノ酸のLC-MS分析

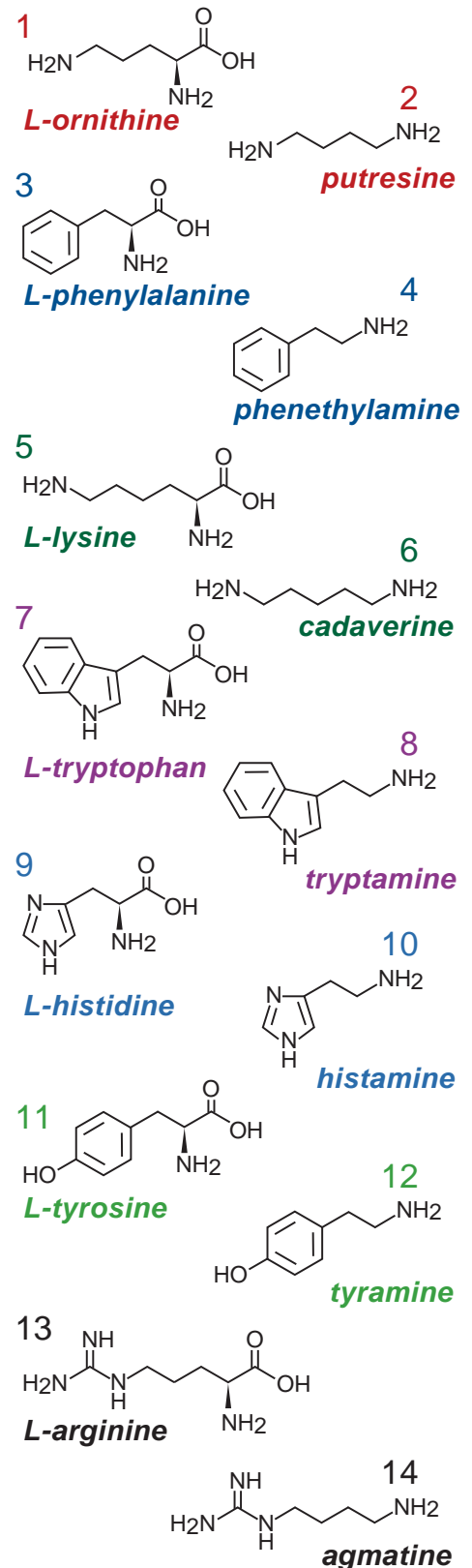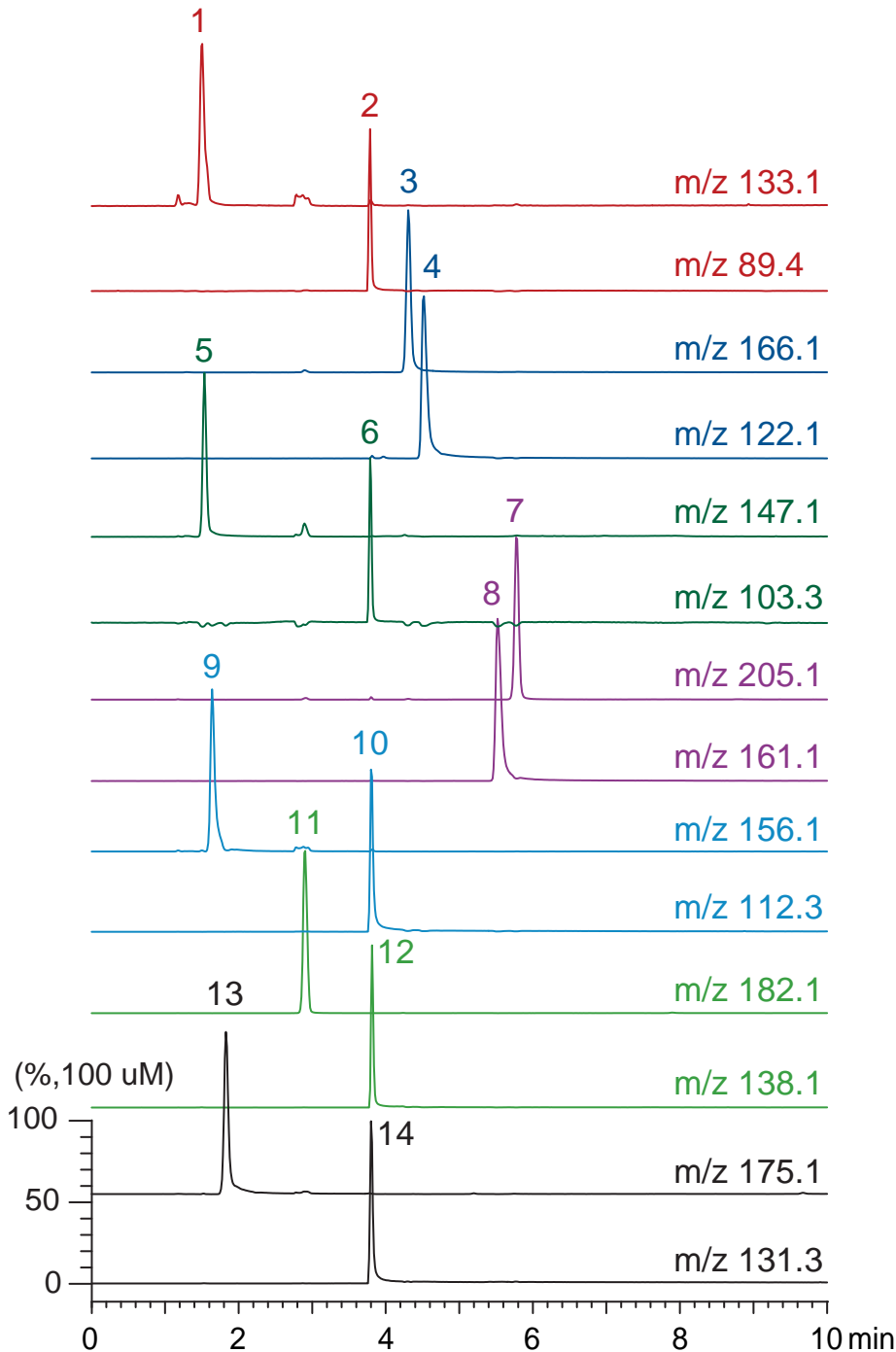

Scherzo SM-C18, 100 x 3 mm

A: 5 mM ammonium formate

B: methanol / water / formic acid = 80 / 20 / 0.5

0-70 %B (0-10min), 100 %B (10-15min)

0.4 mL/min (10 MPa), 37°C, 5uL (100nmol/mL, 0.01N HCl)

Single Quad. MS (ESI, positive)

Distributor

LCC Engineering & Trading GmbH, Steinbruchstr. 4, CH-4622 Egerkingen, Switzerland.

www.chromatographyshop.com , info@chromatographyshop.com,

Tel: +41 62 398 52 71 , Fax: +41 62 398 52 74