



Scherzo SM-C18

150 x 4.6 mm

Application

S-allyl-L-cysteine in Garlic Extract

ニンニク抽出液中のアリルシステイン

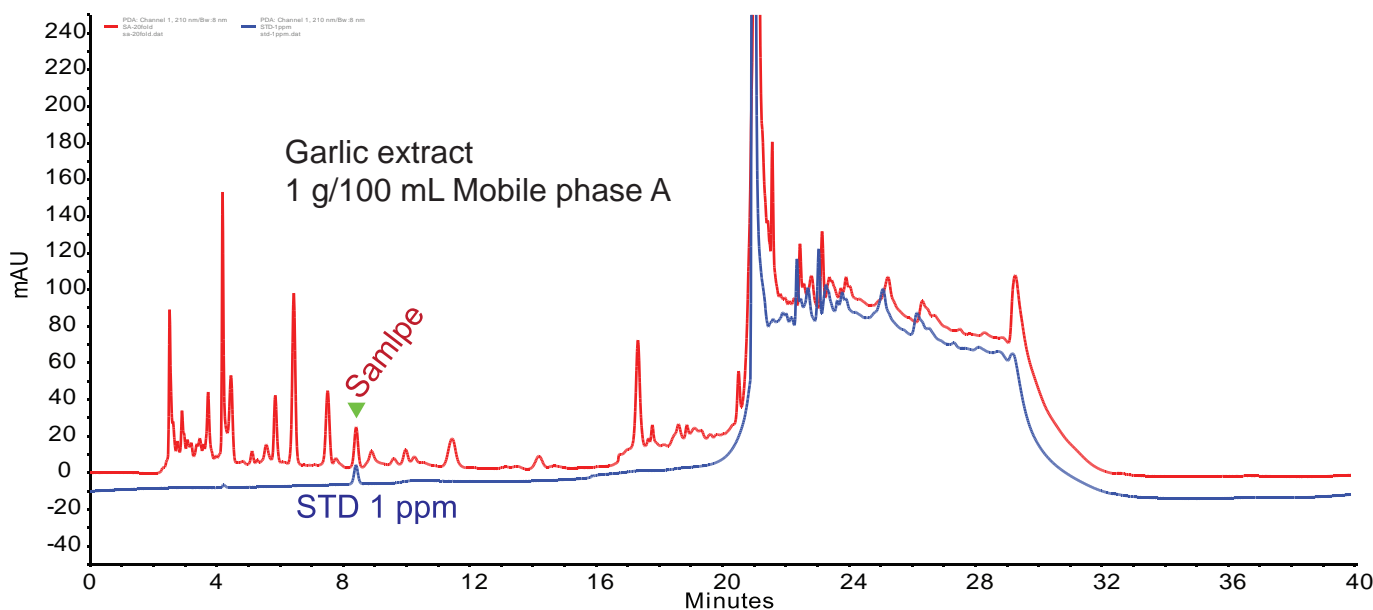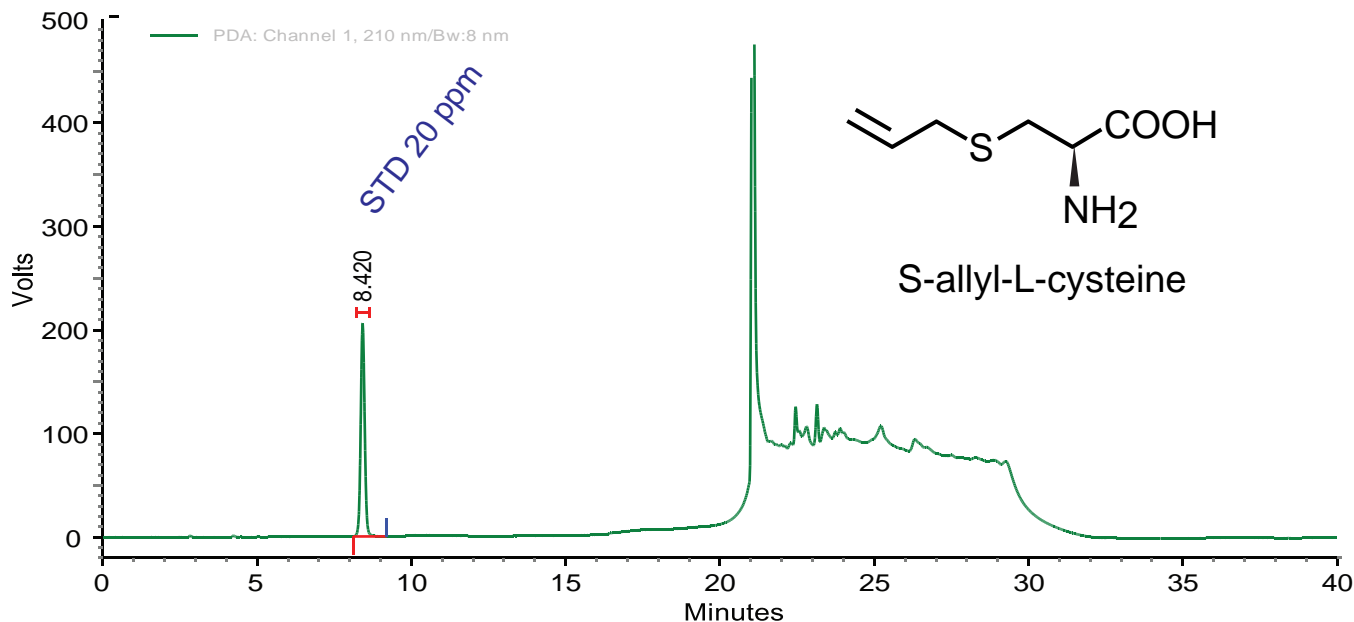

Scherzo SM-C18, 150 x 4.6 mm

A: 0.1% H₃PO₄ in water

B: 0.1% H₃PO₄ in 80% acetonitrile

0%B (0-7min), 0-10%B (7-15min), 10-90%B (15-17min), 90%B (17-25min), 0%B (25-40min)

0.5 mL/min, 40°C

5 µL

210nm

Courtesy of PEAKMAN SP, KOREA

Distributor

LCC Engineering & Trading GmbH, Steinbruchstr. 4, CH-4622 Egerkingen, Switzerland.

www.chromatographyshop.com , info@chromatographyshop.com,

Tel: +41 62 398 52 71 , Fax: +41 62 398 52 74

