

10 Different Cannabinoids and Compounds of Interest

Application #AN6850

Conditions

Column: ACE Excel 2 SuperC18
Dimensions: 100 x 3.0 mm
Part Number: EXL-1011-1003U
Mobile Phase: A: 0.075% formic acid in H₂O
B: 0.1% formic acid in MeCN

| Time (mins) | %B |
|-------------|-----|
| 0.0 | 70 |
| 1.0 | 70 |
| 10.0 | 100 |
| 12.0 | 100 |

Flow Rate: 0.6 mL/min
Injection: 5 µL
Temperature: 40 °C
Detection: UV, 210 nm

1. (-)-11-Nor-9-carboxy- Δ^9 -tetrahydrocannabinol (THC-COOH)
2. Cannabidiarin (CBDV)
3. Cannabidiolic acid (CBDA)
4. Cannabigerol (CBG)
5. Cannabidiol (CBD)
6. Tetrahydrocannabivarin (THCV)
7. Cannabinol (CBN)
8. (-)-*trans*- Δ^9 -Tetrahydrocannabinol (THC)
9. Cannabichromene (CBC)
10. Δ^9 -Tetrahydrocannabinolic acid A (THCA-A)

