


Intrada WP-RP
Presto FF-C18


250 x 4.6 mm

Application

Collagen
コラーゲン

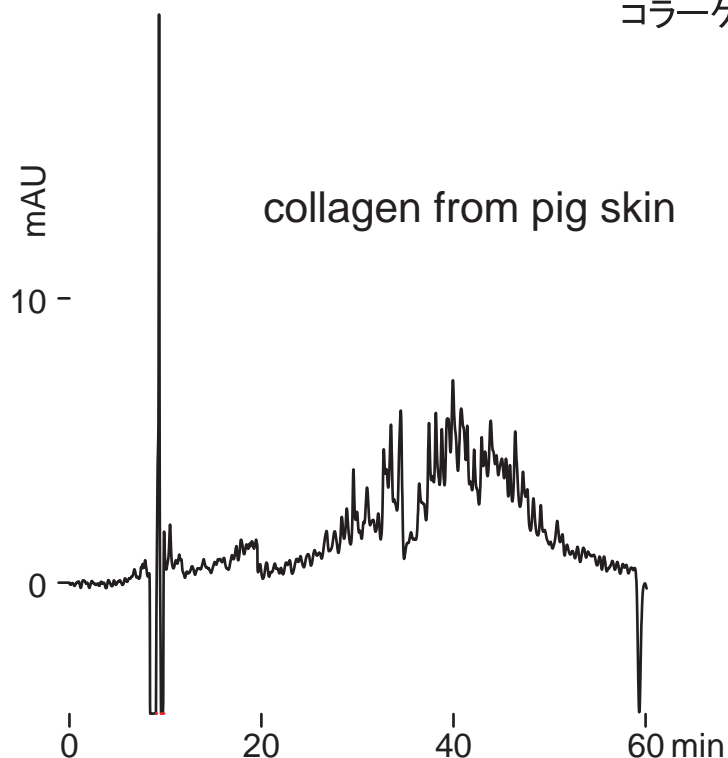

Intrada WP-RP, 250 x 4.6 mm
 A: water /trifluoroacetic acid = 100 /0.1
 B: water /acetonitrile /trifluoroacetic acid = 40 /60 /0.1
 15-25 %B (0-50min)
 0.4 mL/min (3MPa), 80 deg.C, 220 nm
 100 uL (0.01%, 1mM HCl)

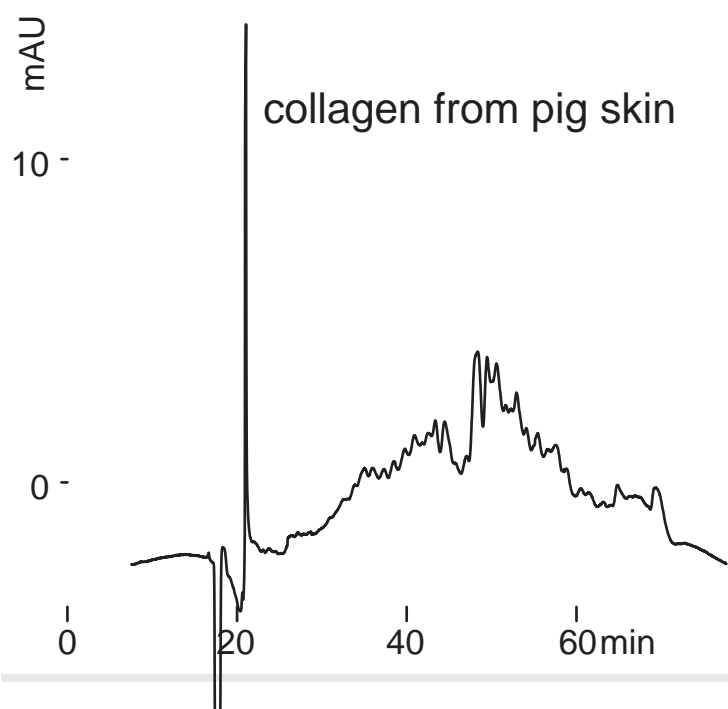

Presto FF-C18, 250 x 4.6 mm
 A: water /trifluoroacetic acid = 100 /0.1
 B: water /acetonitrile /trifluoroacetic acid = 40 /60 /0.1
 35-45 %B (0-50min)
 0.15 mL/min (7MPa), 80 deg.C, 220 nm
 50 uL (0.01%, 1mM HCl)

Available from

LCC Engineering & Trading GmbH
 Steinbruchstrasse 4, CH-4622 Egerkingen, Switzerland
 Tel.: +41 62 398 5271, Fax.: +41 62 398 5274
info@chemsupply.ch
www.chromatographyshop.com

Intakt Corp./JAPAN (info@imtakt.com)
 Other Countries

TI686E-KH29